



**AATCC SERISE**

**AATCC    Dryer machine**

**Model:LBT-M6D**

**Operation instructions**

1/Objective: to use the washing machine properly and correctly

2/Suitable for Laboratory Using

3/Reference standard:

AATCC TM88B、AATCC TM88C、AATCC 124、AATCC 135、AATCC 143、AATCC 150、AATCC LP1-2018

4/Use methods and steps:

1Man-machine interactive console



## 2、Time Setting:




- 1) By manual knob, drying time can be selected according to the demand. Time decreases to the left and increases to the right. Generally, the initial default is 70 minutes drying time.
- 2) Choose the appropriate time according to the demand before drying clothes.


### 3、Drying Procedure Setting:



- 1) The left knob of the panel is to select the drying procedure. There are three procedures: NO Heat (no heating mode), Delicate (low temperature mode) and Normal/Permanent press (high temperature mode).
- 2) The right side of the panel is the display interface of the drying program.

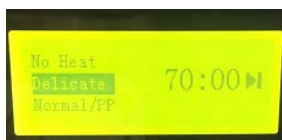

- 3) The drying mode corresponding to the setting of the process knob will be displayed

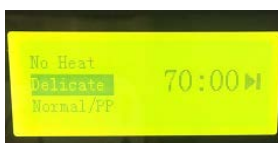


on the display screen.

The drying operation begins:

- 1) When the drying operation begins, the drying mode and time should be set first.



- 2) Put the object to be dried into the barrel and press the key . The dryer will run automatically and the green working indicator will be on.
- 3) When the drying operation is completed, the equipment will stop automatically, and

the red light working instruction will be turned on.

- 4) If you need to stop halfway, you can press the start button and the dryer will stop running.



- 5) If you need to cancel the whole program, press about three seconds, then the operation page of the dryer will return to the main page, and then proceed with the selection of the program.

- 6) The standard program Delicate (low temperature mode), Normal/Permanent press (high temperature mode), has 10 minutes of cold air drying time. When the confirmation program runs, the display page will be added 10 minutes of time to display, while the NO Heat file will be displayed according to the adjusted actual time.